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DE	Gebrauchshinweis – SARSTEDT PCR-Gefäße	2 - 3
EN	Instructions for Use – SARSTEDT PCR vessels	4 - 5
BG	Инструкции за употреба – SARSTEDT PCR епруветки	6 - 7
CS	Návod k obsluze – zkumavky PCR SARSTEDT	8 - 9
DA	Brugsanvisning – SARSTEDT PCR-rør	10 - 11
EL	Οδηγίες χρήσεως – Περιέκτες PCR SARSTEDT	12 - 13
ES	Instrucciones de uso – Microtubos para PCR SARSTEDT	14 - 15
ET	Kasutusjuhend – SARSTEDT PCR-anumad	16 - 17
FR	Mode d'emploi – Tubes PCR SARSTEDT	18 - 19
HR	Uputa za upotrebu – SARSTEDT PCR epruvete	20 - 21
HU	Használati utasítás – SARSTEDT PCR cső	22 - 23
IT	Istruzioni d'uso – Contenitori per PCR SARSTEDT	24 - 25
KO	사용 설명서 – SARSTEDT PCR 튜브	26 - 27
LT	Naudojimo instrukcijos – SARSTEDT PCR mėgintuvėliai	28 - 29
LV	Lietošanas norādes – SARSTEDT PCR trauki	30 - 31
NL	Gebruiksaanwijzing – SARSTEDT PCR-buisjes	32 - 33
NO	Bruksanvisning – SARSTEDT PCR-rør	34 - 35
PL	Instrukcja obsługi – naczynia do PCR SARSTEDT	36 - 37
PT	Instruções de utilização – Tubos SARSTEDT PCR	38 - 39
RO	Instrucțiuni de utilizare – tuburi PCR SARSTEDT	40 - 41
RU	Инструкция по применению – ПЦР-пробирки SARSTEDT	42 - 43
SK	Návod na Použitie – SARSTEDT PCR skúmavky	44 - 45
SL	Navodila za uporabo – epruvete PCR SARSTEDT	46 - 47
SV	Bruksanvisning – SARSTEDT PCR-rör	48 - 49
TH	คำแนะนำในการใช้งาน – หลอดเก็บตัวอย่าง PCR ของ SARSTEDT	50 - 51
TR	Kullanım için talimatlar – SARSTEDT PCR Tüpleri	52 - 53
ZH	使用说明 – SARSTEDT PCR管	54 - 55

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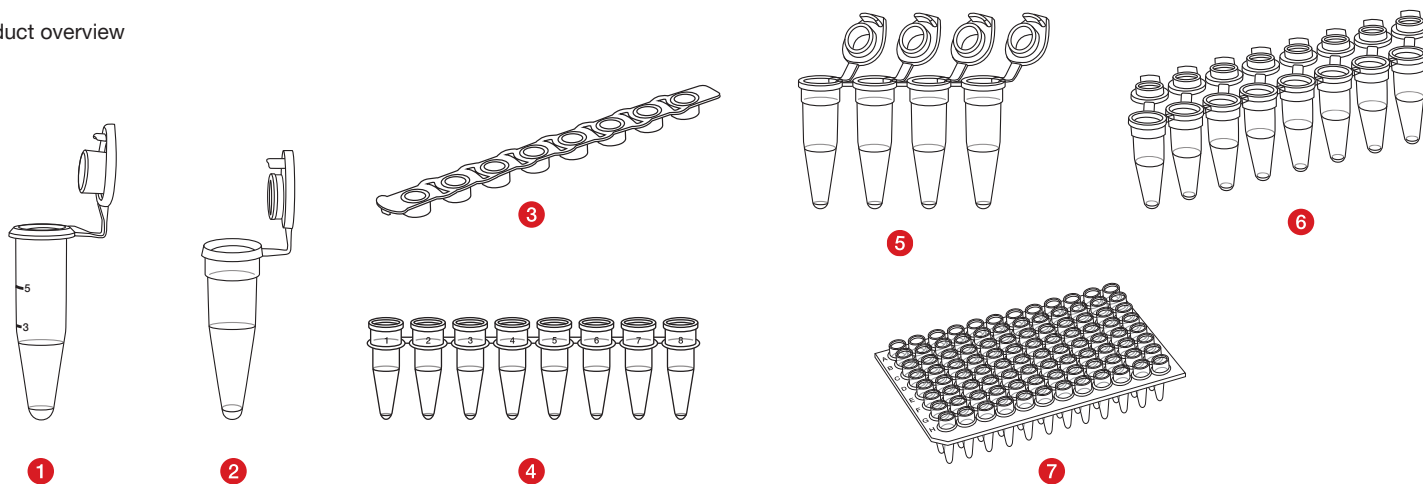
Intended use

PCR vessels are used to hold reaction mixtures, to perform reactions (e.g. PCR) and to store (>24 h) specimen material. PCR vessels are general laboratory supplies. PCR vessels are intended for a single use only by qualified personnel trained in diagnostic PCR applications

Product description

PCR vessels are available in different variants (see product overview). The variants have different colours, shapes, maximum volumes, and cap types. There are also sterile and non-sterile variants and PCR vessels of different purity standards such as PCR Performance Tested and Biosphere® plus, as well as various packaging options available.

Product overview

	Art. no.	Description
1	72.735.xxx	MultiPLY® Pro PCR vessel 0.5 ml
2	72.737.xxx	MultiPLY® Pro PCR vessel 0.2 ml
7	72.985	96-well PCR plate, Biosphere® plus
4	72.985.xxx	MultiPLY® μStrip 0.2 ml, 8 tubes, without cap
5	72.990.xxx	MultiPLY® μStripPro 0.2 ml, 4 tubes
6	72.991.xxx	MultiPLY® μStripPro, 8 tubes

Cap strips:

	Art. no.	Description
3	65.989.xxx	Strip of 8 caps, flat, highly transparent

Safety information and warnings

1. Personal protective equipment must be used when handling diagnostic specimens or hazardous substances.
2. Do not use PCR vessels after the expiry date.
3. Biological specimens must be handled in compliance with the applicable hygiene and safety regulations.
4. The PCR vessels are intended for single use. After use, they must be disposed of in accordance with applicable regulations. There is a risk of contamination if reused.

Storage

The product should be stored at room temperature; avoid exposure to direct sunlight.

Handling

1. Open the bag containing the PCR vessels and take them out.
2. Position container chains for use in a compatible stand.
3. Place your specimen material and any reagents into the PCR vessel. Make sure that the vessel is not filled above the nominal volume.
4. Close the PCR vessel with the attached cap or a separate cap strip.
5. If applicable, transfer the PCR vessels to a thermal cycler and start the PCR program.

Disposal

1. Wearing gloves and other personal protective equipment reduces the risk of infection. Follow the regulations at your facility.
2. The general hygiene guidelines and statutory regulations for the proper disposal of infectious material must be noted and complied with.
3. Contaminated PCR vessels must be disposed of in appropriate biohazard disposal containers. Follow the regulations at your facility.

Key for symbols and labels:

Article number

Batch number

Use by

CE mark

For *In-vitro* diagnostic use

Follow the instructions for use

Risk of contamination if reused

Keep away from sunlight

Do not use if package is damaged

Manufacturer

Country of manufacture

The following is also applicable to sterile products:

Sterilisation by EO

non-pyrogenic

Technical modifications reserved.

All serious incidents relating to the product shall be notified to the manufacturer and the competent national authority.

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