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## SARSTEDT Pipettenspitzen

<b>DE</b>	Gebrauchshinweis – SARSTEDT Pipettenspitzen	2 - 5
<b>EN</b>	Instructions for Use – SARSTEDT Pipette Tips	6 - 9
<b>BG</b>	Инструкции за употреба – SARSTEDT накрайници за пипети	10 - 13
<b>CS</b>	Návod k obsluze – pipetovací špičky SARSTEDT	14 - 17
<b>DA</b>	Brugsanvisning – SARSTEDT pipettespidser	18 - 21
<b>EL</b>	Οδηγίες χρήσεως – Ρύγχη πιπέτας SARSTEDT	22 - 25
<b>ES</b>	Instrucciones de uso – Puntas de pipeta SARSTEDT	26 - 29
<b>ET</b>	Kasutusjuhend – SARSTEDT pipetiotsakud	30 - 33
<b>FI</b>	Käyttöohje – SARSTEDT-pipetinkärjet	34 - 37
<b>FR</b>	Mode d'emploi – Pointes de pipette SARSTEDT	38 - 41
<b>HR</b>	Uputa za upotrebu – SARSTEDT nastavci za pipetu	42 - 45
<b>HU</b>	Használati utasítás – SARSTEDT pipettahegyek	46 - 49
<b>IT</b>	Istruzioni d'uso – Puntali per pipette SARSTEDT	50 - 53
<b>KO</b>	사용 설명서 – SARSTEDT 피펫 팁	54 - 57
<b>LT</b>	Naudojimo instrukcijos – SARSTEDT pipėčių antgaliai	58 - 61
<b>LV</b>	Lietošanas norādes – SARSTEDT pipēšu uzgali	62 - 65
<b>NL</b>	Gebruiksaanwijzing – SARSTEDT-pipetpunten	66 - 69
<b>NO</b>	Bruksanvisning – SARSTEDT pipettespisser	70 - 73
<b>PL</b>	Instrukcja obsługi – końcówki do pipet SARSTEDT	74 - 77
<b>PT</b>	Instruções de utilização – Pontas de pipeta SARSTEDT	78 - 81
<b>RO</b>	Instrucțiuni de utilizare – vârfuli de pipete PCR SARSTEDT	82 - 85
<b>RU</b>	Инструкция по применению – Наконечники для пипеток SARSTEDT	86 - 89
<b>SK</b>	Návod na Použitie – pipetové špičky SARSTEDT	90 - 93
<b>SL</b>	Navodila za uporabo – konice pipet SARSTEDT	94 - 97
<b>SV</b>	Bruksanvisning – SARSTEDT pipettspetsar	98 - 101
<b>TH</b>	คำแนะนำในการใช้งาน – ปิเปตต์ที่ปลายสำหรับหัวของเหลวของ SARSTEDT	102 - 105
<b>TR</b>	Kullanım için talimatlar – SARSTEDT Pipet uçları	106 - 109
<b>ZH</b>	使用说明 – SARSTEDT移液器吸头	110 - 113

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## Intended use

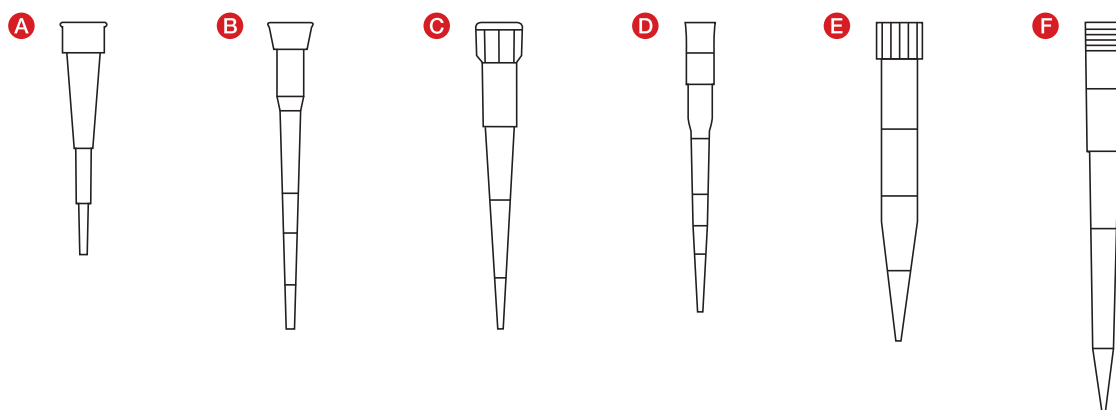
SARSTEDT pipette tips are suitable for a low-contamination transfer of fluids used for *in vitro* diagnostic examinations. The tips are for single use only and intended for use by qualified personnel.

No special ambient conditions are required.

## Product description

The pipettes are available in different variants. The designs vary primarily in pipette volume and their compatibility with the pipettors available on the market that are tested according to DIN EN ISO 8655-6. An overview of the tested combinations (pipette tips – pipettor) can be found at [www.sarstedt.com](http://www.sarstedt.com). In addition, different packaging variants, pipette tips with or without filter, tips with a low retention surface, and many different purity standards such as PCR Performance Tested and Biosphere® plus are available.

## Product overview

## Pipette tips:

Type	Art. no.	Description
<b>A</b>	70.3010.xxx	10 µl pipette tip
<b>B</b>	70.3020.xxx / 70.3021.xxx	20 µl pipette tip
<b>C</b>	70.3030.xxx / 70.3031.xxx	200 µl pipette tip
<b>D</b>	70.3040.xxx	300 µl pipette tip
<b>E</b>	70.3050.xxx	1000 µl pipette tip
<b>F</b>	70.3060.xxx	1250 µl pipette tip

## Safety information and warnings

1. Personal protective equipment must be used when handling diagnostic specimens or hazardous substances.
2. Do not use pipette tips after the expiry date.
3. Biological specimens must be handled in compliance with the applicable hygiene and safety regulations.
4. The pipette tips are intended for single use only. After use, they must be disposed of in accordance with applicable regulations. There is a risk of contamination if reused.

## Storage

Store the pipette tips at room temperature and avoid exposure to direct sunlight.

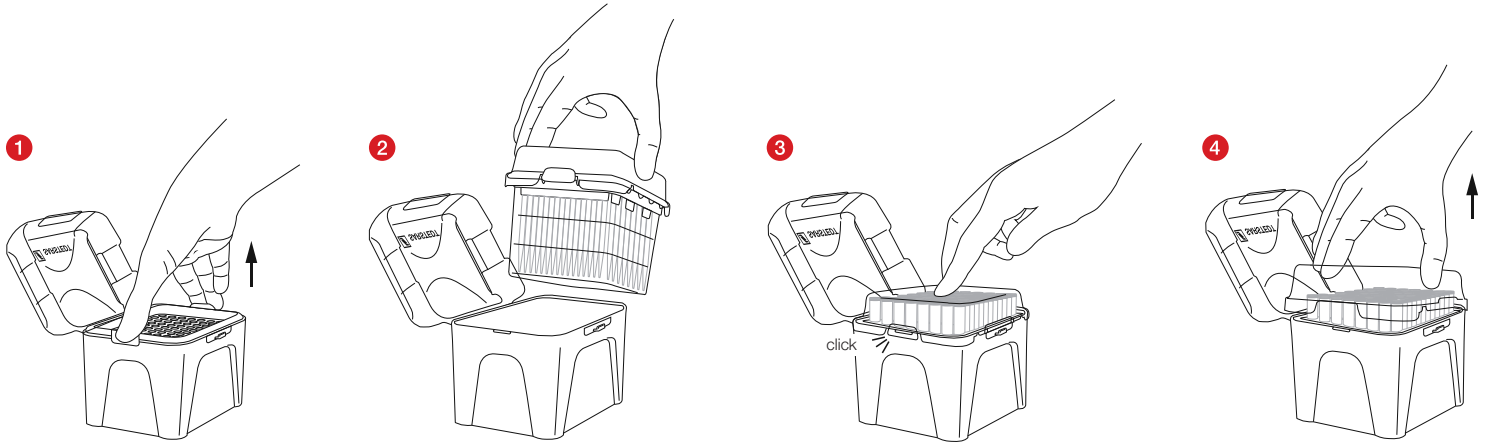
## Limitations

For use only with pipettors that have been tested for compatibility. An overview of the tested combinations (pipette tips – pipettor) can be found at [www.sarstedt.com](http://www.sarstedt.com).

**Handling**

1. Open the packaging of the pipette tip box and take it out.
2. Open the lid of the box by pressing lightly on "PRESS".
3. Attach the pipette tip to your pipettor.


**Refilling the box with SingleRefill:**

- 1 Release the empty tray by pressing on the clips at the sides simultaneously.
- 2 Transfer the SingleRefill to the box.
- 3 Press on the middle of the lid: the SingleRefill clicks into place with an audible "CLICK".
- 4 Now simply lift up the lid to remove it.

**Refilling the box with StackPack:**

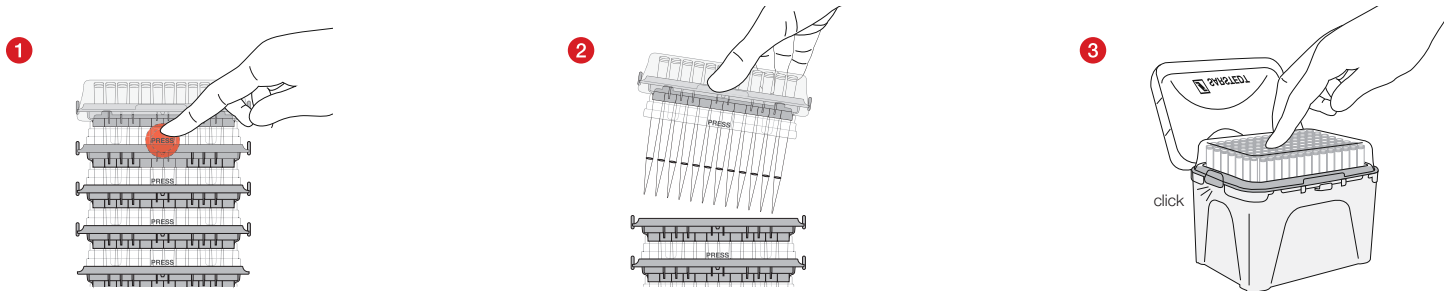
**Variant 1**


- 1 Press the "PRESS" points at the front and back one after the other.
- 2 Lift up the single tray to remove it and transfer it to the box.
- 3 Press lightly on the middle of the lid until it clicks firmly into place.

Variant 2

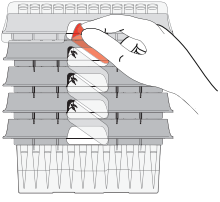1

2

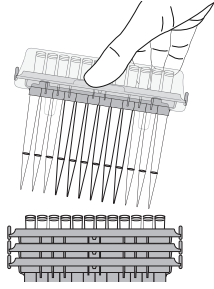3

- 1 Remove the uppermost label on each side one after the other.
- 2 Lift up the single tray to remove it and transfer it to the box.
- 3 Press lightly on the middle of the lid until it clicks firmly into place.

Disposal

1. Wearing gloves and other personal protective equipment reduces the risk of infection. Follow the regulations at your facility.
2. The general hygiene guidelines and statutory regulations for the proper disposal of infectious material must be noted and complied with.
3. Contaminated pipette tips must be disposed of in appropriate biohazard disposal containers. Follow the regulations at your facility.

US standards / ISO standards (as amended / updated)

Product-specific standards

DIN EN ISO 8655-2: Piston-operated volumetric apparatus - Part 2: Piston pipettes

DIN EN ISO 8655-6: Piston-operated volumetric apparatus - Part 6: Gravimetric reference test method for determining volume

Key for symbols and labels:

	Article number
	Batch number
	Use by
	CE mark
	For <i>in-vitro</i> diagnostic use
	Follow the instructions for use
	Risk of contamination if reused
	Keep away from sunlight
	Store in a dry place
	Manufacturer
	Country of manufacture

The following is also applicable to sterile products:

	ETO sterile
	Single sterile barrier system with external protective packaging

Technical modifications reserved.

All serious incidents relating to the product shall be notified to the manufacturer and the competent national authority.

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